

PCR Check System

Quantity: 20 reactions

Store: -20°C (not frost free)

Cat. No.: PK3091

Description: The *PCR Check system* provides qualified reagents for the control of PCR reaction. PCR Check System is supplied with quality tested Taq DNA Polymerase (#DP1602), 10X PCR buffer (#MM2101), and a vial of MgCl₂ (#MM2091), 10mM dNTP mix (#MM2082), DNA template and related "Reverse" and "Forward" Primers. For added convenience, nucleotides are provided as a 10mM dNTP mix. The *PCR Check system's* reagents are sufficient for 20 reactions of 25µl each.

Components (supplied):

Taq DNA Polymerase (5 unit/µl)	8µl
10X PCR Buffer	50µl
50mM Magnesium Chloride	15µl
dNTP Mix (10mM)	10µl
Control DNA (1pg/µl)	100µl
Forward Primer	20µl
Reverse primer	20µl
Distilled Water	5ml
Mineral Oil	1ml

Control DNA: (Supplied) DNA extracted from an insert in plasmid. This DNA has been thoroughly tested for its performance in several PCR reactions with various annealing temperatures ranging from 40-65°C. The amplified DNA shows a band of 620bp in an Agarose gel. This DNA may be used in parallel as control.

Control forward and Reverse primer:

(Supplied) These primers have successfully been used and tested in several PCR reactions with mentioned control DNA.

PCR Runs: The *PCR Check system* is used for amplification of different reactions.

Data obtained from all of the reactions was satis-factory.

Basic PCR protocol:

The following basic protocol serves as a general guideline and a starting point for any PCR amplification. Optimal reaction conditions (incubation times, temperatures, concentration of Taq DNA Polymerase (#DP1602), primers, MgCl₂, and template DNA) vary and to be evaluated by the user. If you have any problem with your PCR system, you can check each reagent by *PCR Check System* kit and evaluate your PCR system.

Add the following components to a sterile 0.5ml micro-centrifuge tube sitting on ice:

Components.....Volume.....Final Con.

Autoclaved distilled water	14µl	
10X PCR buffer	2.5µl	0.2mM each
10mM dNTP mix	0.5µl	0.2mM each
50mM MgCl ₂	0.75µl	1.5mM
Forward primer	1µl	0.5µM
Reverse primer	1µl	0.5µM
Control DNA (4pg/µl) 5µl	20pg	
Taq DNA polymerase	0.2µl	1 unit
Mineral Oil	25µl	

Steps Temperature Range of duration (seconds)

Denature	95°C		120sec.	
Denature	93°C	10sec.	45sec.	} 35 cycles
Anneal	40-65°C	10sec.	60sec.	
Extend	72°C	20sec.	90sec.	

Result analysis:

Analyze 10µl of amplified samples in a 2% agarose gel after adding loading buffer. The presence of 620bp fragment indicates of positive.

Centrifuge Tubes Before Opening To Improve Recovery Of Contents.

For Research Use Only

SinaClon BioScience

www.sinaclon.com

Tel: +98(0)21 4463 0050-51

+98(0)21 4466 5156

Customer Service: +98(0)21 4463 3016

Place your order: order@sinaclon.com